
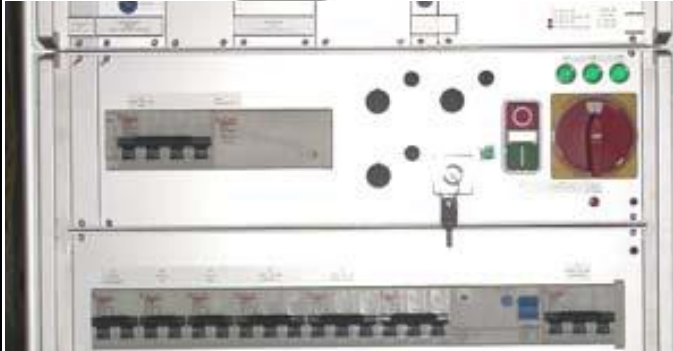


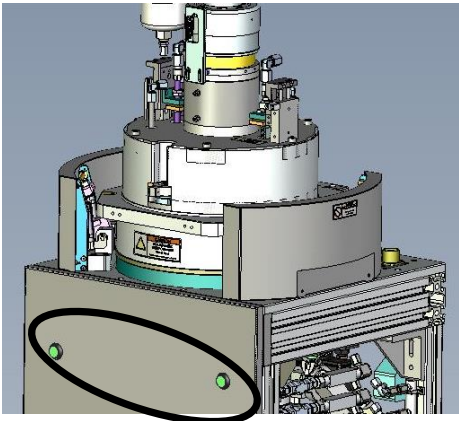


## OPERATING INSTRUCTIONS EVG 520 HOT EMBOSSE

	<p>Turn On MAIN SWITCH (RED) of the embosser by rotating in clockwise direction</p> <p>The three green light on the top should come on</p> <p>Push the GREEN BUTTON on the switch left of MAIN SWITCH</p> <p>The PC has no separate switch and should turn on now</p>
	<p>Push the two GREEN BUTTONS at the front of the process chamber to open/close the chamber and load your sample</p> <p>Note: The buttons should be pressed until the chamber is either fully open or fully closed. In fully closed mode, the clips surrounding the chambers will come down and lock the chamber. Fully open means until it does not open further. You can also read the status of the lid in the software. If the position is in between, it will say UNDEFINED</p>
<ul style="list-style-type: none"><li>• Login by clicking button at the bottom left</li><li>• Click on  button at the bottom to edit a recipe</li><li>• You can change the embossing parameters such as time, temperature, pressure etc and SAVE the recipe</li><li>• Click on  to LOAD and RUN the recipe</li><li>• Click on  . It will ask you to select the recipe and then prompt for wafer and lot ID</li></ul>	

- Once the process is finished, it will display message that the process has finished. Click on LOAD NEXT to load another sample with same recipe or END to exit.

### **SHUTTING DOWN**

- Exit the program by clicking “FILE” -> “EXIT”
- Shutdown Windows
- Press the RED button of switch next to the MAIN SWITCH
- Turn OFF MAIN SWITCH