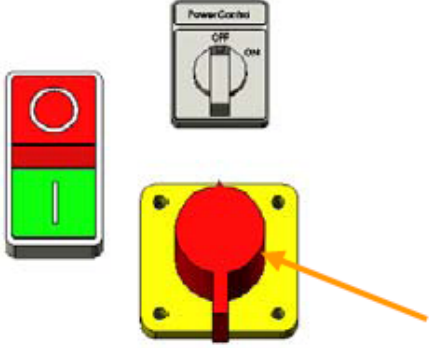

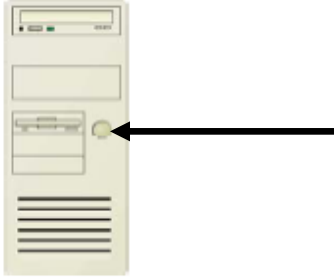


# OPERATING INSTRUCTIONS EVG 620 MASK-ALIGNER

	<p>Turn On MAIN SWITCH of the mask aligner by rotating in clockwise direction</p> <p>Push the GREEN BUTTON on the switch left of MAIN SWITCH</p> <p>Note: The KEY-SWITCH of the electronic must be OFF; otherwise you can't ignite the lamp</p>
	<p>Turn ON POWER SWITCH of lamp power supply</p> <p>Press START SWITCH on the lamp power supply (left hand side)</p> <p>Wait about 10 minutes until the lamp has reached the working temperature and the whole power before starting the exposure. However, you can start the electronics immediately after the lamp has ignited.</p>
	<p>Turn KEY-SWITCH to ON position to power the electronics</p>
	<p>Turn ON PC SWITCH of the PC and login</p>
<ul style="list-style-type: none"> <li>• Press "FILE" -&gt; "OPEN" to open a saved recipe</li> <li>• Edit the exposure parameters such as TYPE, TIME etc and save the recipe by "FILE" -&gt; "SAVE"</li> <li>• Click on RUN to start the recipe and follow the instructions on the screen</li> </ul>	

- After all the exposures are finished involving the same mask, unload the mask by clicking on UNLOAD MASK button on the recipe window. Follow the instructions on the screen
- Once the mask is unloaded, click on PARK TRAY to move the tray in and follow the instructions on the screen.

### **SHUTTING DOWN**

- Exit the program by clicking “FILE” -> “EXIT”
- Shutdown Windows
- Turn OFF Electronic KEY-SWITCH
- Turn OFF Lamp Power Supply
- Press POWER SWITCH of the Lamp Power Supply unit for cooling
- Wait 10 minutes
- Turn OFF Lamp Power Supply
- Press the RED button of switch next to the MAIN SWITCH
- Turn OFF MAIN SWITCH